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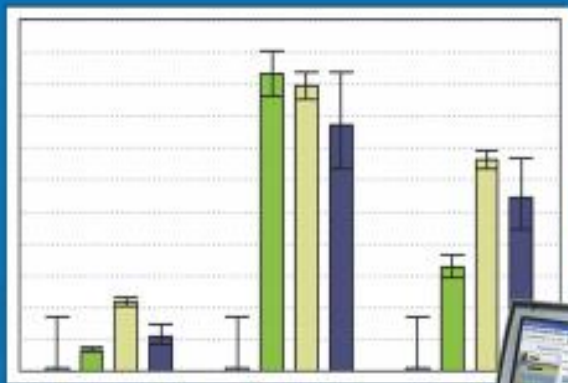
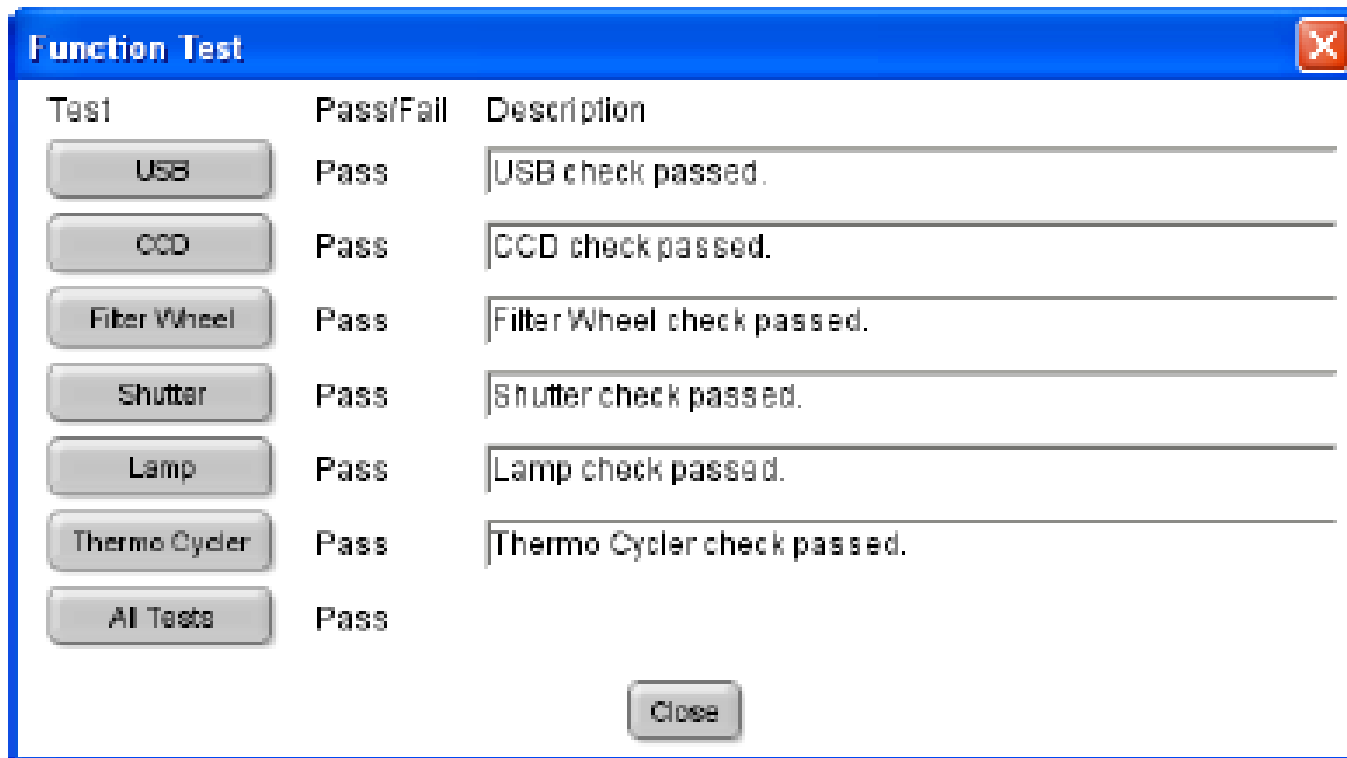
The screenshot shows a software window titled "Function Test" with a close button in the top right corner. The window contains a table with three columns: "Test", "Pass/Fail", and "Description". Each row represents a different component that has been tested and found to pass. At the bottom of the window, there is a "Close" button.


Test	Pass/Fail	Description
USB	Pass	USB check passed.
CCD	Pass	CCD check passed.
Fiber Wheel	Pass	Filter Wheel check passed.
Shutter	Pass	Shutter check passed.
Lamp	Pass	Lamp check passed.
Thermo Cycler	Pass	Thermo Cycler check passed.
All Tests	Pass	

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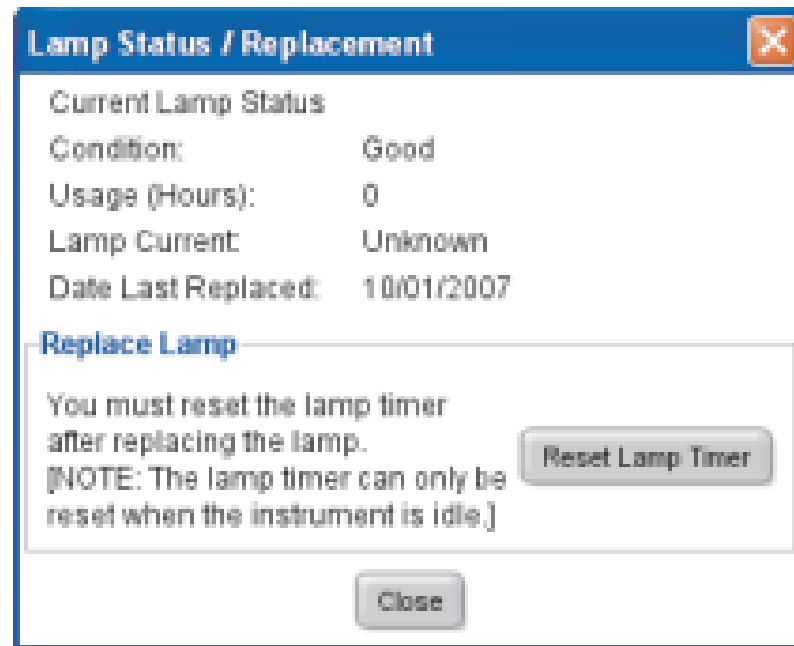**Lamp Status / Replacement** 

Current Lamp Status

Condition:	Good
Usage (Hours):	0
Lamp Current:	Unknown
Date Last Replaced:	10/01/2007

**Replace Lamp**

You must reset the lamp timer after replacing the lamp.  
[NOTE: The lamp timer can only be reset when the instrument is idle.]

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
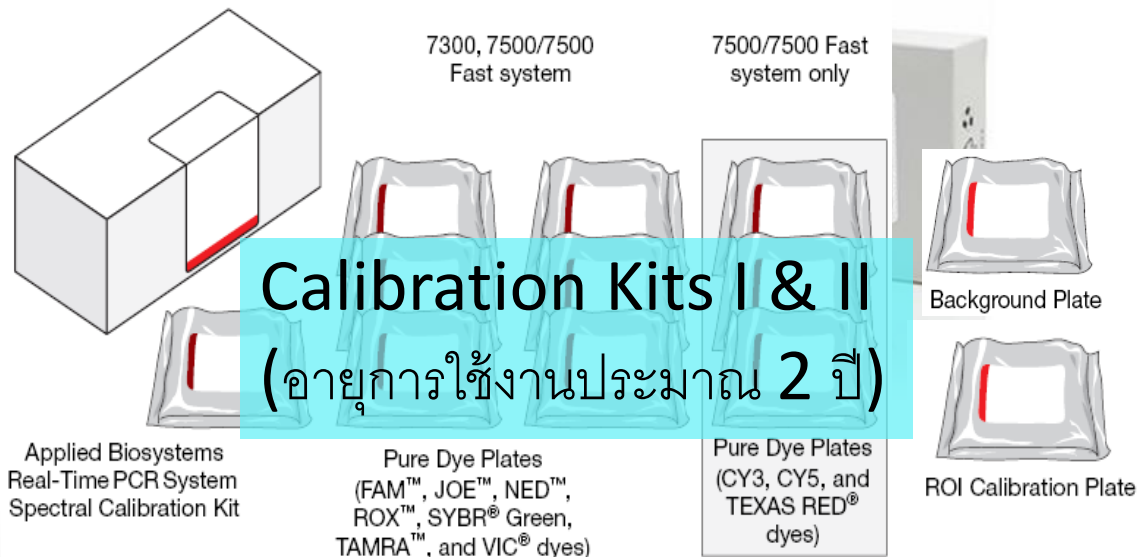
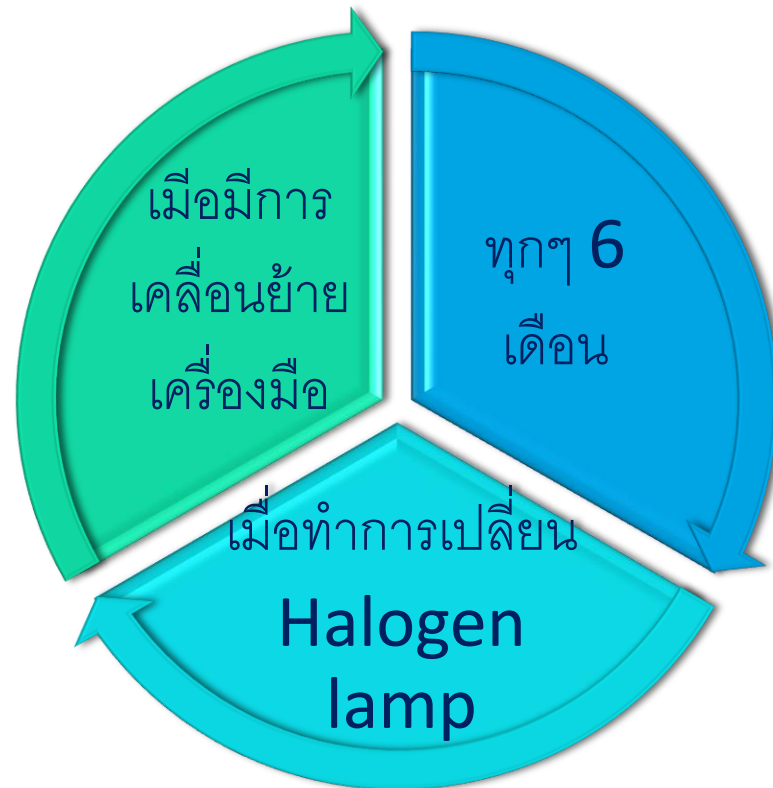
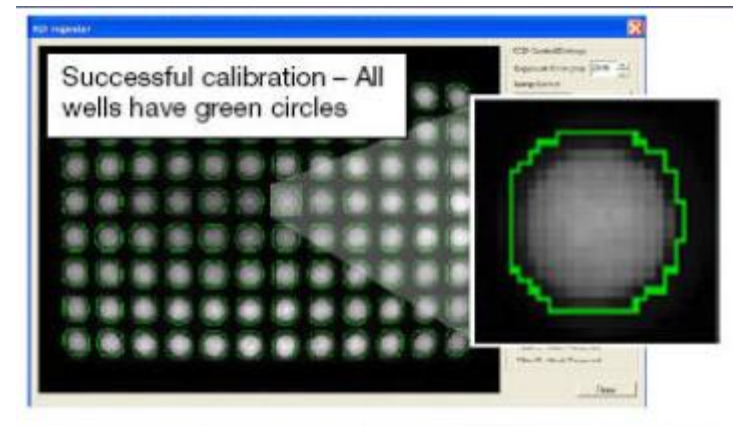
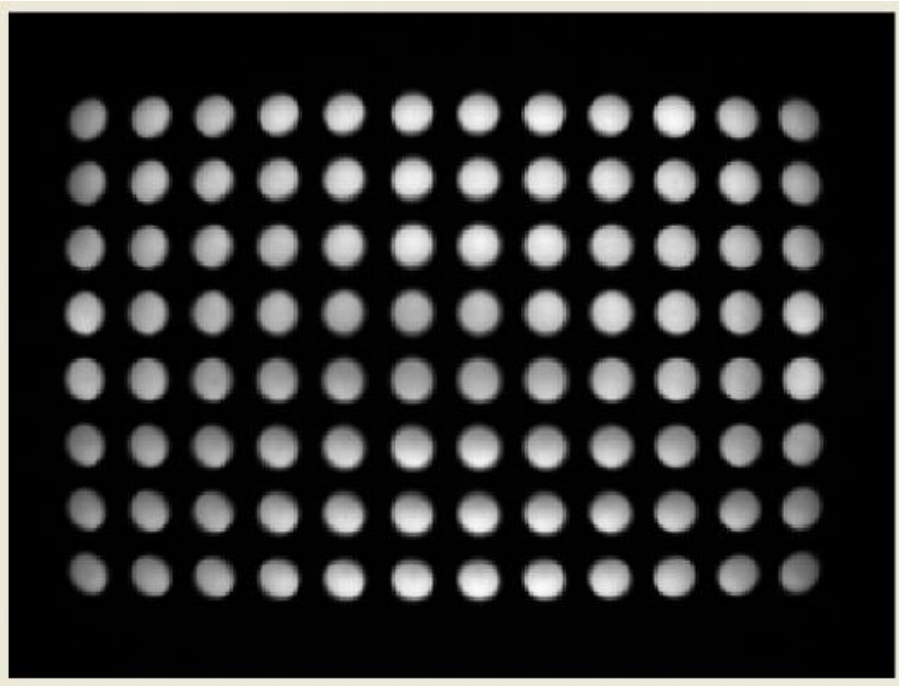
# Region of Interest (ROI) calibration

Load standard plates (7500 system) with the notched A12 position at the top-right of the tray.

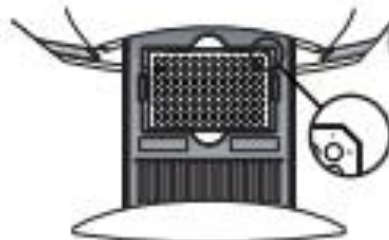Load Fast plates (7500 Fast system) with the notched A1 position at the top-left of the tray.

Close the tray door. Apply pressure to the right side of the tray door at an angle.

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

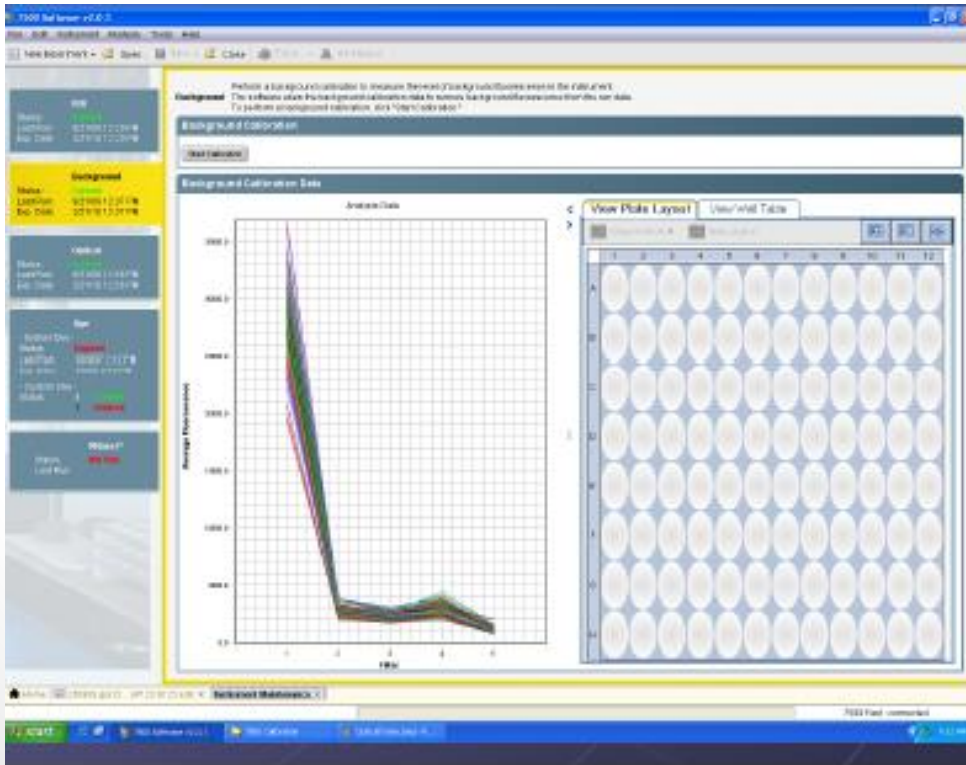
# Sample block decontamination

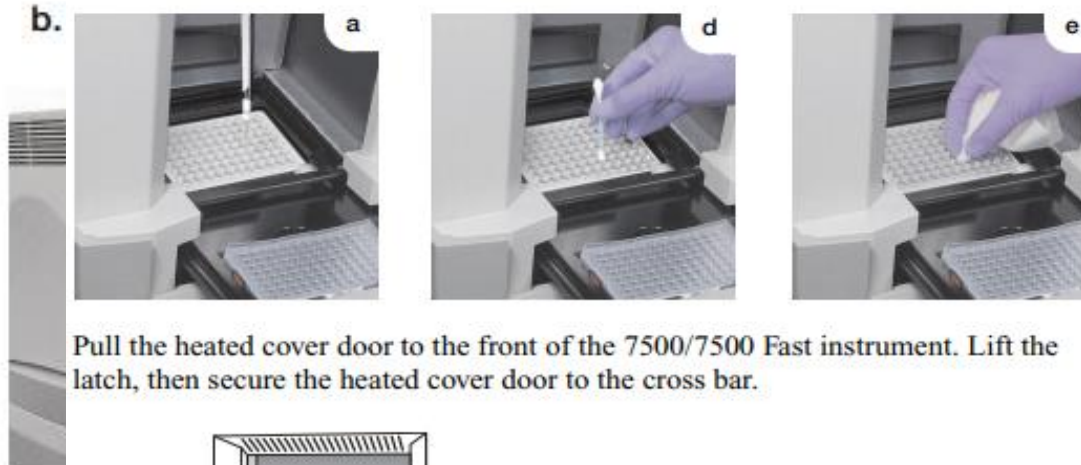- d. Using a cotton swab, scrub inside of each contaminated well.
- e. Using a lint-free cloth, absorb the excess deionized water.

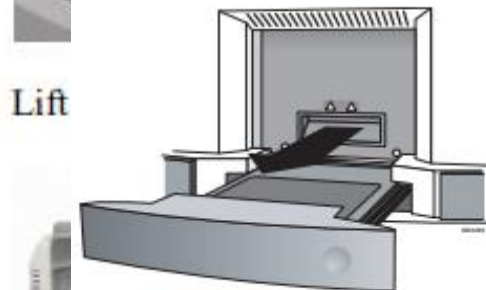Pull the heated cover door to the front of the 7500/7500 Fast instrument. Lift the latch, then secure the heated cover door to the cross bar.

Lift


Close the access door to the 7500/7500 Fast inst

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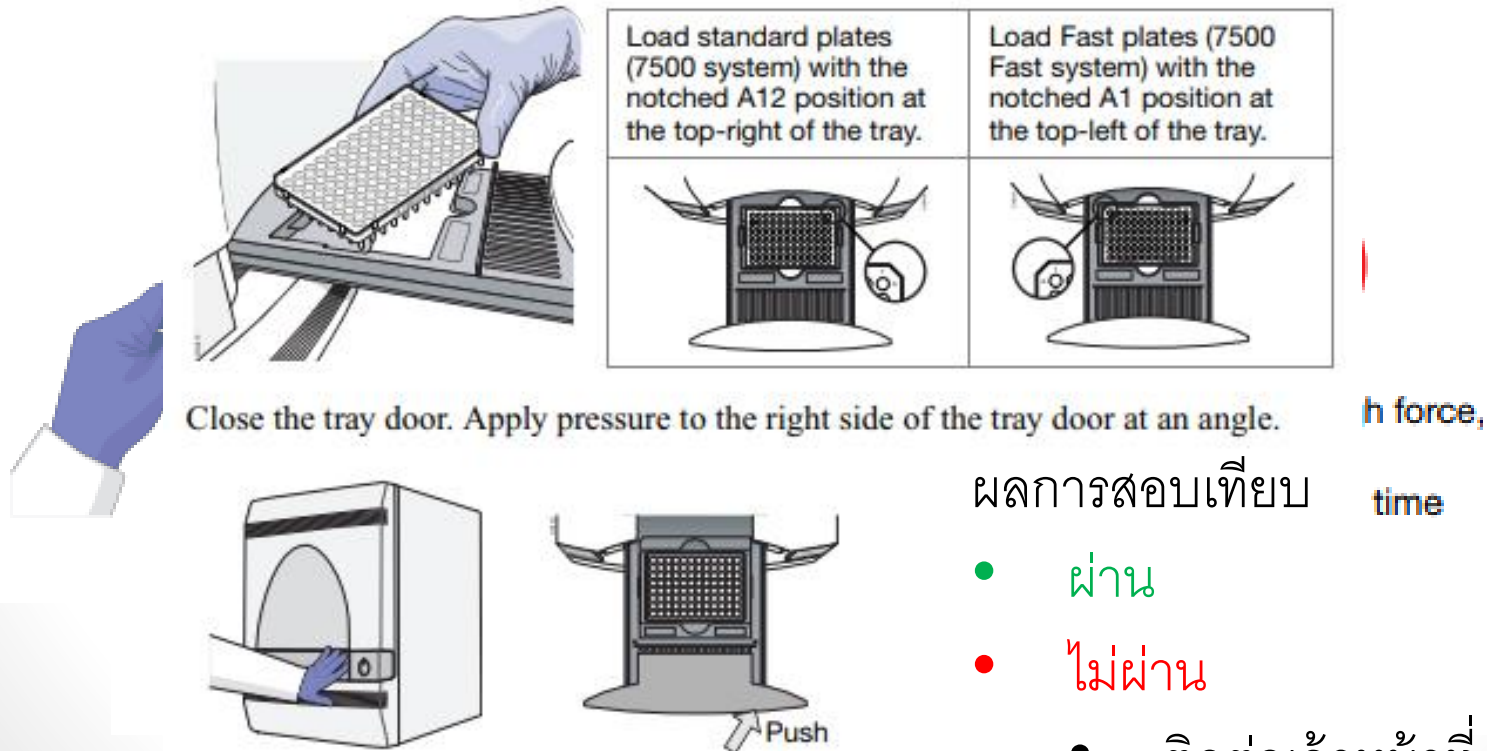
# Pure Dyes calibration

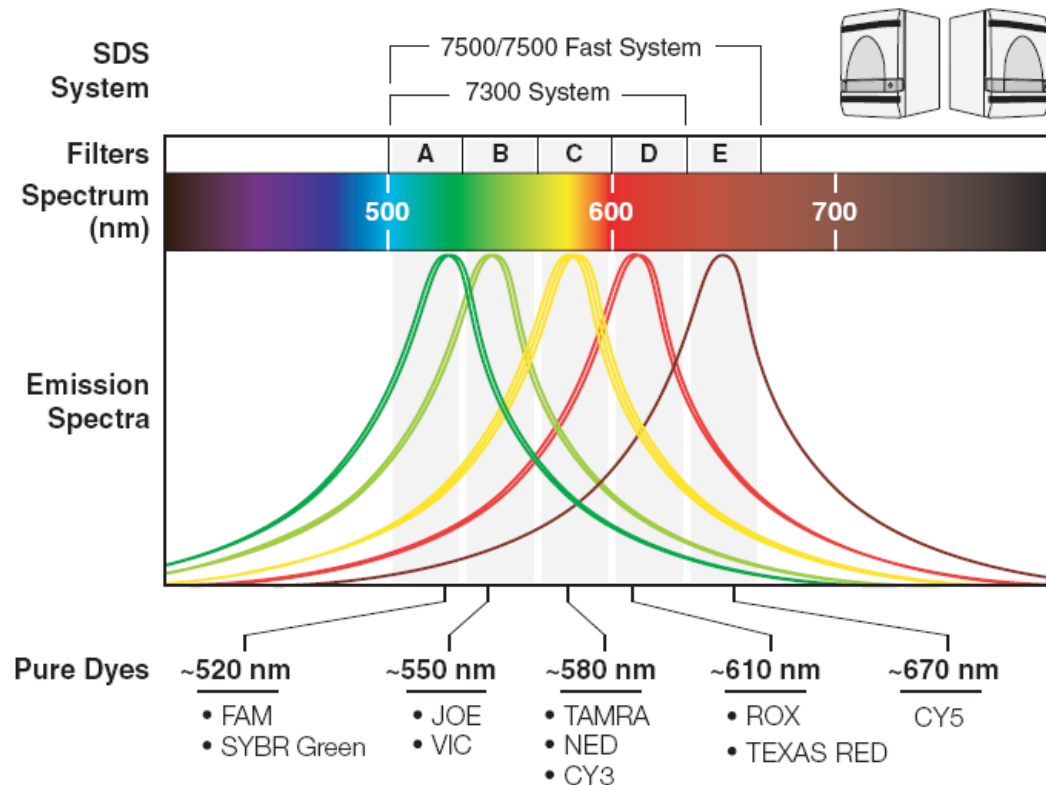- Spectral Information for each of the pure dyes is stored
- Allows the instrument to determine the individual contribution of each dye from the fluorescence collected during a run containing multiple dyes

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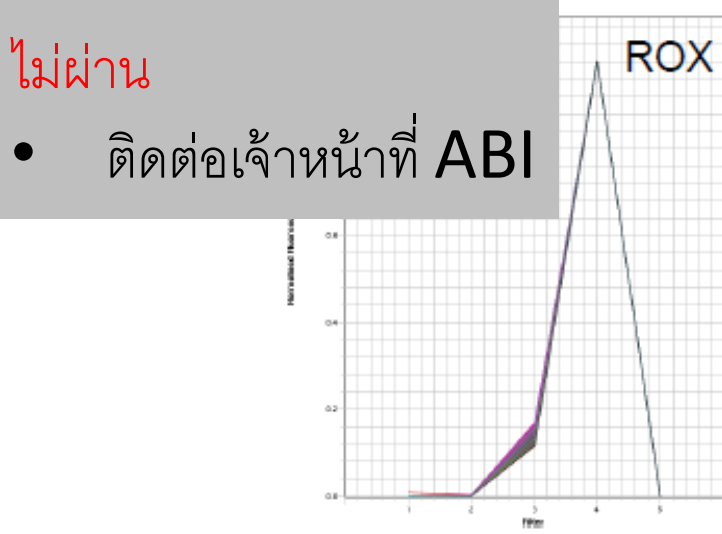
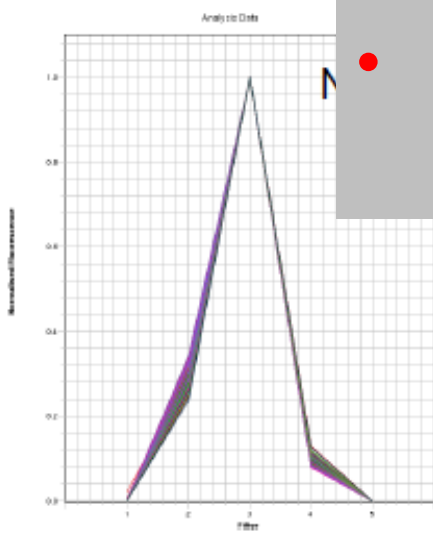
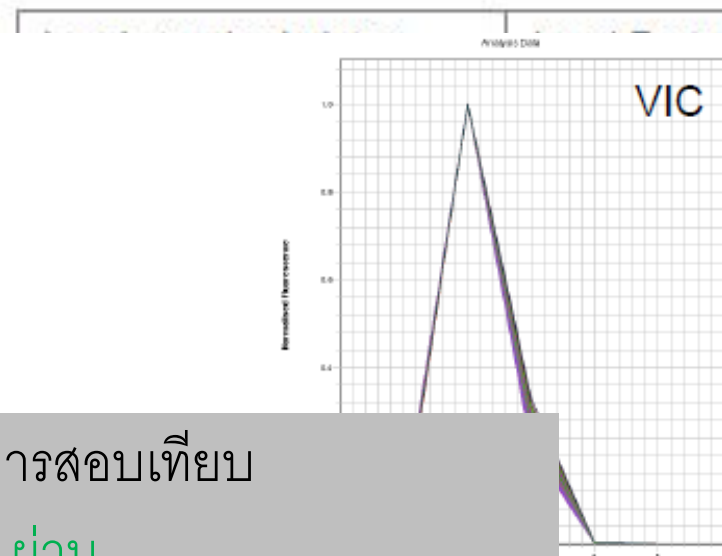
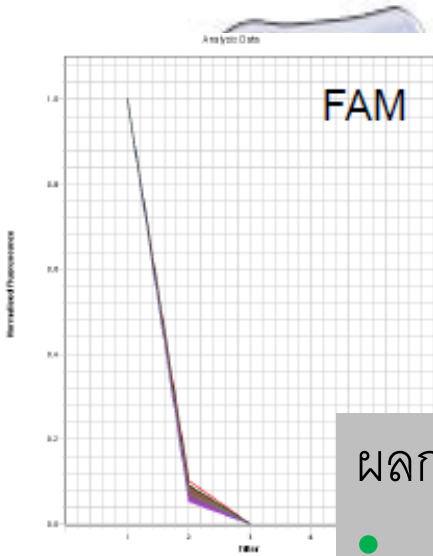
# RNase P calibration

**RNase P Run** Analysis Help

**Instructions:** Review the results from the RNase P run. If the run failed, omit outliers, then click "Reanalyze." Repeat until the run passes. Click **Finish** to save and apply the results.

Status: Passed

Buttons: Omit Reanalyze Analysis Settings

Amplification Plot Standard Curve

Plot Settings

Target: All Plot Color: Default

Save current settings as the default

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